



thermo scientific

QuEChERS solutions in all your favorite flavors



Find it at eu.fishersci.com/go/thermochrom

 **fisher scientific**
part of Thermo Fisher Scientific

QuEChERS: Simple, proven, sample extraction and clean-up

Fast, easy and cost-effective, QuEChERS is a robust, reproducible approach for extracting and cleaning up pesticides and other low-level contaminants in complex matrices. It is often used as a sample preparation step prior to GC-MS or LC-MS analysis, and can be automated for higher throughput.

Originally developed in 2003 for the determination of multiple pesticide residues in food, QuEChERS applications have evolved to include:

- Pesticide residues
- Veterinary drug residues
- Pharmaceutical residues
- Environmental contaminants
- Mycotoxins
- Drugs of abuse
- Traditional Chinese medicines
- Forensic analysis




Pre-packaged, ready-weighed kits make QuEChERS easier and more convenient

Thermo Scientific™ QuEChERS kits give you reproducible results and excellent recoveries for a wide variety of analytes, and save time and money, too. Pre-packaged, ready-weighed salts, solid-phase extraction (SPE) sorbents and buffers streamline your workflow, and minimize the potential for error.

Kits are available in several formats to meet all your application requirements. Depending on your analytes of interest, your sample matrix and your preferred method—Original, AOAC or EN—you can choose the appropriate volume option, and select the right combination of salts, sorbents and consumables.



QuEChERS workflow

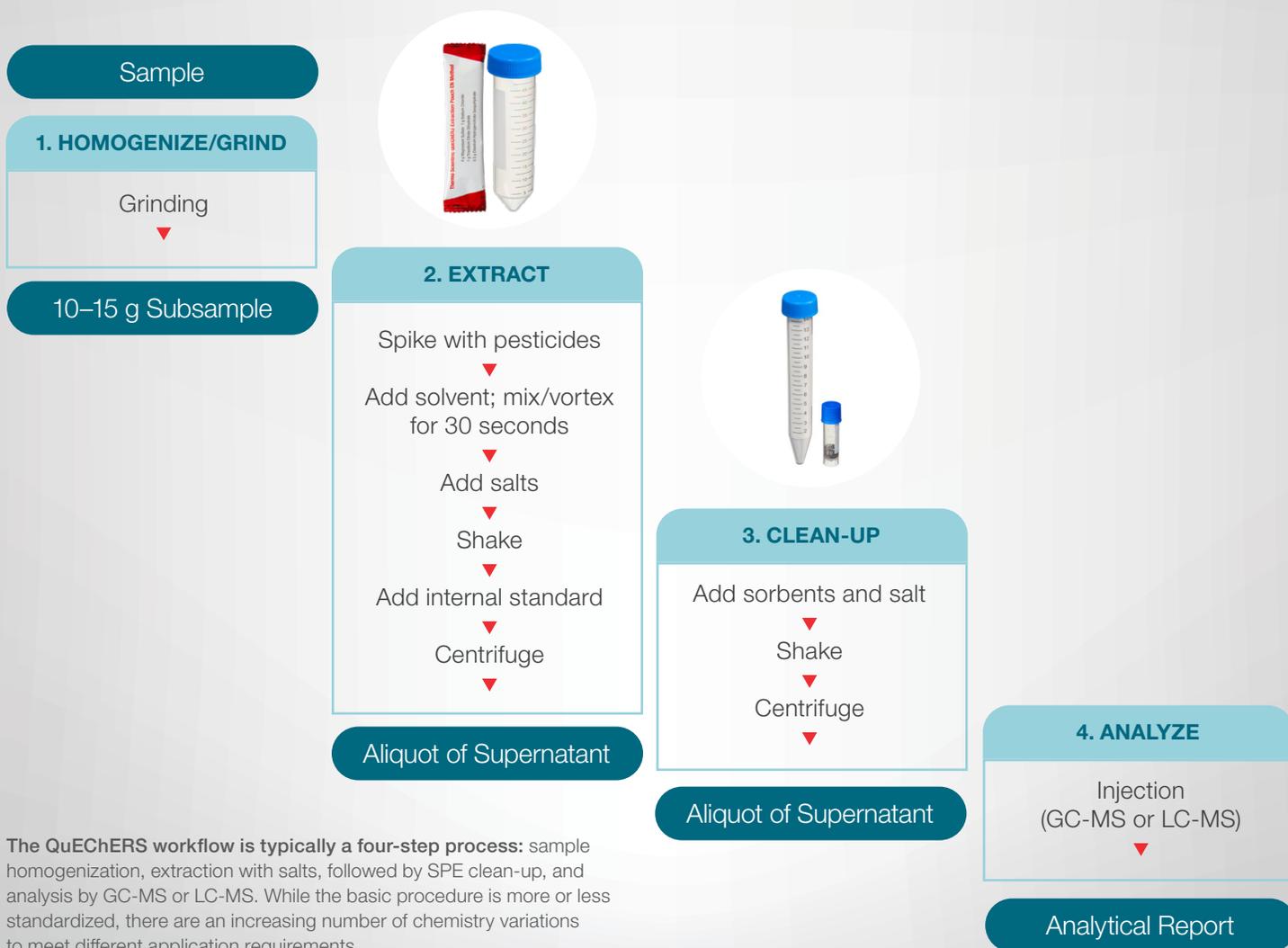

The QuEChERS workflow is typically a four-step process: sample homogenization, extraction with salts, followed by SPE clean-up, and analysis by GC-MS or LC-MS. While the basic procedure is more or less standardized, there are an increasing number of chemistry variations to meet different application requirements.

Three flavors of QuEChERS

Kits and products to support them all

Whichever QuEChERS method you're using, we've got you covered

Original

(Anastassiades et al, 2003):

developed for non-base-sensitive compounds, using sodium chloride to reduce polar interferences and enhance extraction

AOAC

(AOAC 2007.01):

uses sodium acetate as a buffer instead of sodium chloride and is compatible with base-sensitive compounds

European

(EN 15662):

similar to the AOAC method, but uses sodium chloride to minimize polar interferences, and sodium citrate dihydrate and disodium citrate sesquihydrate instead of sodium acetate



Thermo Scientific QuEChERS Extraction Pouch EN Method

4 g Magnesium Sulfate 1 g Sodium Chloride
1 g Trisodium Citrate Dihydrate
0.5 g Disodium Hydrogencitrate Sesquihydrate



Original method kits

Original method extraction kits*

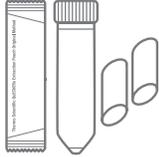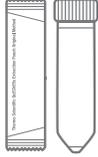Extract	Description	Capacity	Quantity	Cat. No
	 <p>Original extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouch): 4 g MgSO₄, 1 g NaCl • 50 mL tubes (empty) • Ceramic homogenizers 	10 g samples	50/PK	17399822
	 <p>Original extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouch): 4 g MgSO₄, 1 g NaCl • 50 mL tubes (empty) 	10 g samples	50/PK	17309832
	 <p>Original extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouch): 4 g MgSO₄, 1 g NaCl 	10 g samples	50/PK	17319832


*Original method extraction kits are also available in 15 g sample capacity.

Bulk ceramic homogenizers — all methods

Homogenize/Grind	Description	Quantity	Cat. No
	 <p>Ceramic homogenizers for 2 mL tubes</p>	100/PK	17339822
	 <p>Ceramic homogenizers for 15 mL tubes</p>	100/PK	17329822
	 <p>Ceramic homogenizers for 50 mL tubes</p>	100/PK	17349822

AOAC method kits

AOAC method extraction kits


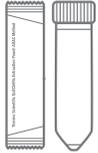
Extract	Description	Capacity	Qty.	Cat. No
	 <p>AOAC extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 6 g MgSO₄, 1.5 g NaCl • 50 mL tubes (empty) • Ceramic homogenizers 	15 g samples	50/PK	17339832
	 <p>AOAC extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 6 g MgSO₄, 1.5 g NaCl • 50 mL tubes (empty) 	15 g samples	50/PK	17349832
	 <p>AOAC extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 6 g MgSO₄, 1.5 g NaCl 	15 g samples	50/PK	17359832

AOAC method clean-up kits

Clean-up	Description	Qty.	Cat. No	
	General fruits and vegetables			
	AOAC Clean-up kit, prefilled 2 mL tubes with 50 mg PSA, 150 mg MgSO ₄	100/PK	17389852	
	AOAC Clean-up kit, prefilled 15 mL tubes with 400 mg PSA, 1200 mg MgSO ₄	50/PK	17329842	
	Pigmented fruits and vegetables			
	AOAC Clean-up kit, prefilled 2 mL tubes with 50 mg PSA, 50 mg GCB, 150 mg MgSO ₄	100/PK	17329862	
	AOAC Clean-up kit, prefilled 15 mL tubes with 400 mg PSA, 400 mg C18, 1200 mg MgSO ₄	50/PK	17329852	
	Fruits and vegetables with fats and waxes			
	AOAC Clean-up kit, prefilled 2 mL tubes with 50 mg PSA, 50 mg C18, 150 mg MgSO ₄	100/PK	17369852	
	AOAC Clean-up kit, prefilled 15 mL tubes with 400 mg PSA, 400 mg C18, 1200 mg MgSO ₄	50/PK	17399832	
	Fruits and vegetables with pigments and fats			
	AOAC Clean-up kit, prefilled 2 mL tubes with 50 mg PSA, 50 mg C18, 50 mg GCB, 150 mg MgSO ₄	100/PK	17349862	
	AOAC Clean-up kit, prefilled 15 mL tubes with 400 mg PSA, 400 mg C18, 400 mg GCB, 1200 mg MgSO ₄	50/PK	17349852	
	All food types			
	AOAC Clean-up kit, prefilled 2 mL tubes with 50 mg PSA, 50 mg C18, 7.5 mg GCB, 150 mg MgSO ₄	100/PK	17359852	
	AOAC Clean-up kit, prefilled 15 mL tubes with 400 mg PSA, 400 mg C18, 45 mg GCB, 1200 mg MgSO ₄	50/PK	17369832	
	Other foods			
	AOAC Clean-up kit, prefilled 2 mL tubes with 25 mg C18, 150 mg MgSO ₄	100/PK	17319862	
	AOAC Clean-up kit, prefilled 15 mL tubes with 150 mg C18, 900 mg MgSO ₄	50/PK	17399842	

EN method kits

EN method extraction kits

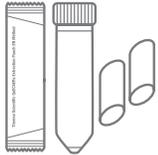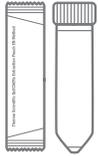	Description	Capacity	Qty.	Cat. No
Extract	 <p>EN extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 4 g MgSO₄, 1 g NaCl, 0.5 g disodium hydrogencitrate sesquihydrate, 1 g trisodium citrate dihydrate • 50 mL tubes (empty) • Ceramic homogenizers 	10 g samples	50/PK	17369822
	 <p>EN extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 4 g MgSO₄, 1 g NaCl, 0.5 g disodium hydrogencitrate sesquihydrate, 1 g trisodium citrate dihydrate • 50 mL tubes (empty) 	10 g samples	50/PK	17379822
	 <p>EN extraction kit</p> <ul style="list-style-type: none"> • Extraction salts (in pouches): 4 g MgSO₄, 1 g NaCl, 0.5 g disodium hydrogencitrate sesquihydrate, 1 g trisodium citrate dihydrate 	10 g samples	50/PK	17389822

EN method clean-up kits


	Description	Qty.	Cat. No
Clean-up	General fruits and vegetables		
	EN Clean-up kit, prefilled 2 mL tubes with 150 mg MgSO ₄ , 25 mg PSA	100/PK	17399852
	EN Clean-up kit, prefilled 15 mL tubes with 900 mg MgSO ₄ , 150 mg PSA	50/PK	17339842
	Pigmented fruits and vegetables		
	EN Clean-up kit, prefilled 2 mL tubes with 150 mg MgSO ₄ , 25 mg PSA, 2.5 mg GCB	100/PK	17339862
	EN Clean-up kit, prefilled 15 mL tubes with 900 mg MgSO ₄ , 150 mg PSA, 15 mg GCB	50/PK	17339852
	Highly pigmented fruits and vegetables		
	EN Clean-up kit, prefilled 2 mL tubes with 150 mg MgSO ₄ , 25 mg PSA, 7.5 mg GCB	100/PK	17309862
	EN Clean-up kit, prefilled 15 mL tubes with 900 mg MgSO ₄ , 150 mg PSA, 45 mg GCB	50/PK	17379842
	Fruits and vegetables with fats and waxes		
EN Clean-up kit, prefilled 2 mL tubes with 150 mg MgSO ₄ , 25 mg PSA, 25 mg C18	100/PK	17379852	
EN Clean-up kit, prefilled 15 mL tubes with 900 mg MgSO ₄ , 150 mg PSA, 150 mg C18	50/PK	17309842	

What is the purpose of the chemicals?

- **MgSO₄**: removes residual water, and induces phase separation between water content in sample and acetonitrile layer
- **NaCl**: removes residual water, and induces phase separation between water content in sample and acetonitrile layer
- **NaOAc**: buffers the sample to stabilize pH
- **Disodium hydrogencitrate sesquihydrate**: buffers the sample to stabilize pH
- **Trisodium citrate dihydrate**: buffers the sample to stabilize pH
- **PSA**: removes free fatty acids and other acidic co-extractives
- **C18**: removes fats, sterols, and other non-polar interferences from sample
- **GCB**: removes pigment (not recommended for use with planar pesticides)

Thermo Scientific VetDrugs Explorer


High-confidence quantitation workflow for veterinary drug residues

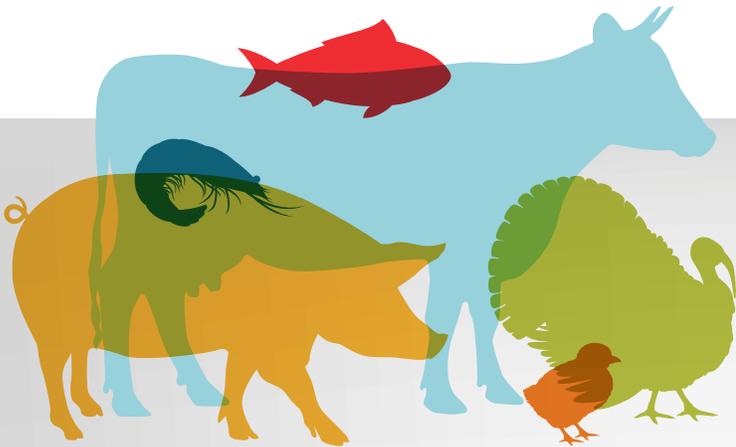



Designed to meet the needs of food safety laboratories aiming for the highest confidence and ultimate sensitivity, the **Thermo Scientific™ VetDrugs Explorer collection** is a validated multi-residue veterinary drug analysis solution that addresses global regulatory requirements.

The end-to-end analytical workflow includes:

- Quality control (QC) and reference standard mixes
- QuEChERS kit with sample preparation procedure
- Industry-leading triple quadrupole mass spectrometer with state of the art UHPLC and columns
- Data acquisition and processing methods, with an extensive compound database—tested and validated by industry thought leaders in multiple laboratories

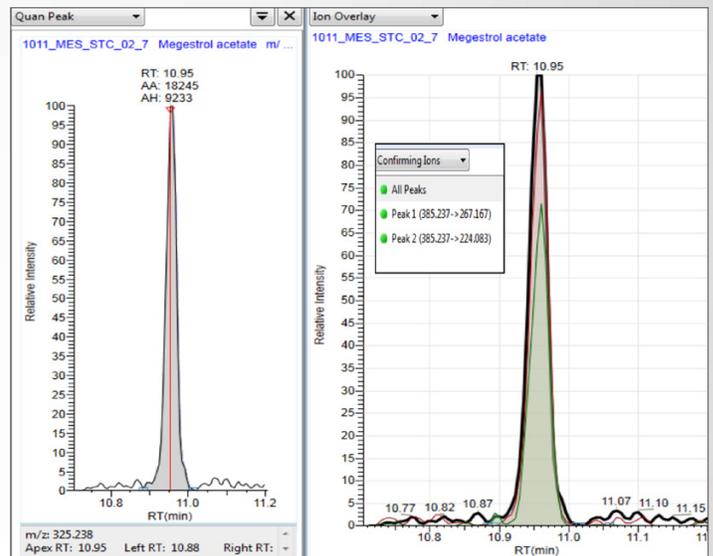

Confident quantitation. Steroid hormone megestrol acetate in milk sample at 0.04 ng/g, quantitation peak (left) with confirming ions (right).

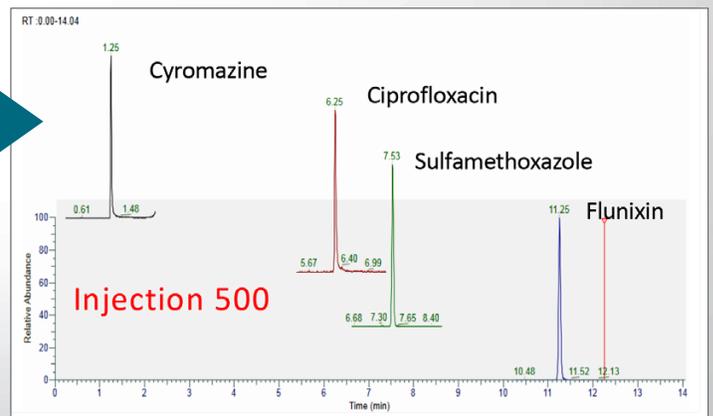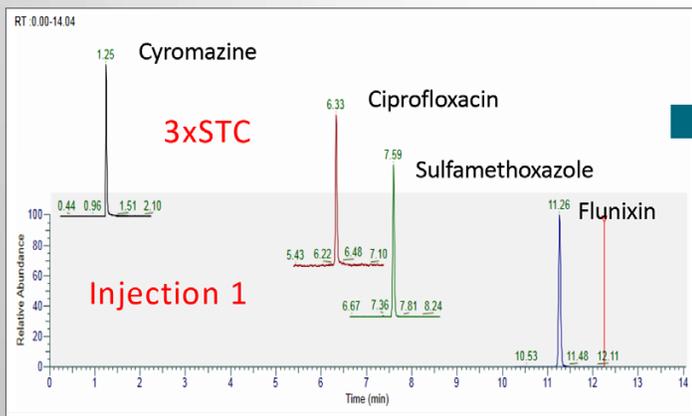

High throughput robustness. Stable peak shape and response over 500 injections in bovine muscle extract.

Today's most comprehensive portfolio of analytical instrumentation

Thermo Scientific™ TSQ series triple quadrupole mass spectrometers

Best-in-class sensitivity and speed, plus outstanding robustness and reliability provide fast, accurate quantitation of hundreds of analytes in a variety of matrices.



Thermo Scientific™ TSQ Altis™ triple quadrupole mass spectrometer



Thermo Scientific™ Orbitrap™ Exploris™ GC 240 mass spectrometer

Thermo Scientific™ Orbitrap™ mass analyzers for GC-MS, LC-MS and MS/MS

High-resolution accurate-mass (HRAM) data allows confident trace-level screening and quantitation of pesticides and identification of unknowns.



Thermo Scientific™ Vanquish™ Horizon UHPLC system

Thermo Scientific™ Vanquish™ UHPLC systems

Today's food safety scientists need a robust system that can support their regulatory analyses. The Vanquish systems improve performance and repeatability with no trade-offs in quality, robustness, or ease-of-use. This innovative, simple-to-operate, and easy-to-maintain platform delivers confident separations of complex food matrices.



Thermo Scientific™ Chromeleon™ Chromatography Data System (CDS)

Proven, industry-leading software streamlines workflows from chromatography to routine quantitative MS analysis on IC, GC, LC, GC-MS/MS and LC-MS/MS systems by delivering superior instrument control, automation, data processing, and more.

Thermo Scientific™ Chromeleon™ Chromatography Data System (CDS)

μSPE Clean-up solution for Thermo Scientific™ TriPlus™ RSH SMART autosampler

Micro solid-phase extraction (μSPE) clean-up solution for the TriPlus RSH SMART autosampler offers a fully automated on-line approach for clean-up of QuEChERS extracts, and replaces the dispersive SPE manual procedure. The TriPlus RSH SMART autosampler is available as a standalone instrument or can be fully integrated with GC-MS or LC-MS systems.



Thermo Scientific™ TriPlus™ RSH SMART autosampler

Recommended consumables

Description	Quantity	Cat. No
GC Method		
QuEChERS GC μSPE Cartridge for use with TriPlus RSH SMART autosampler μSPE, QuEChERS Blend for GC, 45 mg (MgSO ₄ , PSA, C18EC, Carbon)	108/PK	60101-45GC
2 mL clear Snap-It glass vial, 11 mm wide opening	100/PK	17324033
11 mm AS Snap-It Seal Star pre-slit caps	100/PK	17334023
LC Method		
QuEChERS LC μSPE Cartridge for use with TriPlus RSH SMART autosampler μSPE, QuEChERS Blend for LC, 30 mg (Z-Sep, C18, CarbonX)	108/PK	60101-30LC
2 mL clear Snap-It glass vial, 11 mm wide opening	100/PK	17324033
11 mm AS Snap-It Seal Star pre-slit caps	100/PK	17334023

Everything else you'll need for QuEChERS success

In addition to convenient, ready-to-run QuEChERS kits, you'll find all the other ingredients for productive QuEChERS analysis—including pipettors and pipette tips, syringe filters, and industry standard Thermo Scientific™ Sorvall™ benchtop centrifuges. And, of course, a full selection of lab-proven chromatography consumables.



Chromatography consumables

Product Description	Method	Application	Cat. No
Columns			
Thermo Scientific™ LinerGold™: Siltek six baffle PTV liner	GC*	Pesticide residues	10300875
Thermo Scientific™ LinerGold™: SSL splitless liner, single taper, deactivated	GC*	Pesticide residues	15447990
Thermo Scientific™ TraceGOLD™ TG-5SIL MS (30 m x 0.25 mm i.d. x 0.25 µm)	GC*	Pesticide residues	10177894
Thermo Scientific™ Accucore™ aQ C18 Polar Endcapped (100 x 2.1 mm, 2.6 µm)	LC	Pesticide residues	11387481
Thermo Scientific™ Acclaim™ Trinity P1 (anionic pesticides) (100 x 3.0 mm, 3.0 µm)	LC	Pesticide residues	11371943
Thermo Scientific™ Accucore™ RP-MS (100 x 2.1 mm, 2.6 µm)	LC	Vet drug residues	11337531
Thermo Scientific™ Accucore™ aQ C18 Polar Endcapped (100 x 2.1 mm, 2.6 µm)	LC	Mycotoxins	11387481
Vials and caps			
2 mL clear screw thread glass vial	GC/LC	All	17343893
2 mL clear screw thread PP vial (recommended for polar pesticides)	GC/LC	All	17303923
2 mL amber screw thread glass vial	GC/LC	All	17373893
Screw cap for 2 mL screw thread vial. White silicone/red PTFE septa	GC/LC	All	17303873
Solvents and reagents			
Thermo Scientific™ UHPLC-MS grade acetonitrile	GC/LC	All	15329865
Thermo Scientific™ UHPLC-MS grade methanol	GC/LC	All	15319865
Thermo Scientific™ UHPLC-MS grade water	GC/LC	All	15339865
Fisher Chemical™ Optima™ LC-MS grade formic acid	GC/LC	All	10596814
Fisher Chemical™ Optima™ Acetonitrile for HPLC and GC	GC/LC	All	10061044
Fisher Chemical™ Optima Methylene Chloride for HPLC and GC-MS	GC/LC	All	10500341
Fisher Chemical™ Optima 95% n-Hexane for HPLC and GC-MS	GC/LC	All	16504344
Fisher Chemical™ Certified AR, Sodium sulfate anhydrous	GC/LC	All	10746292

*For the complete list of GC consumables for these applications, please refer to our GC Consumables Selection Guide

Find out more at eu.fishersci.com/go/thermochrom

Distributed by Fisher Scientific. Contact us today:

Austria: fishersci.at Belgium: fishersci.be Denmark: fishersci.dk
 Germany: fishersci.de Ireland: fishersci.ie Italy: fishersci.it
 Finland: fishersci.fi France: fishersci.fr Netherlands: fishersci.nl
 Norway: fishersci.no Portugal: fishersci.pt Spain: fishersci.es
 Sweden: fishersci.se Switzerland: fishersci.ch UK: fishersci.co.uk

© 2021 Thermo Fisher Scientific Inc. All rights reserved.

Trademarks used are owned as indicated at fishersci.com/trademarks. BR000249-FEU

